

# CASA<sup>®</sup>

(Campylobacter Selective Agar)

Selective Chromogenic Agar  
for Detection and Enumeration  
of *Campylobacter* spp


## Direct enumeration of *Campylobacter* on CASA<sup>®</sup> plates

- High selectivity: No interfering flora
- Easy reading: brick red colonies
- Simple and easy confirmation with a latex agglutination test

## Detection of *Campylobacter* according to ISO 10272

- All reagents needed for campylobacter detection are available:
  - Bolton broth
  - mCCDA medium
  - CASA<sup>®</sup> medium\*

\*In complement of mCCDA medium, CASA medium provides high selectivity and easy reading, in compliance with ISO 10272

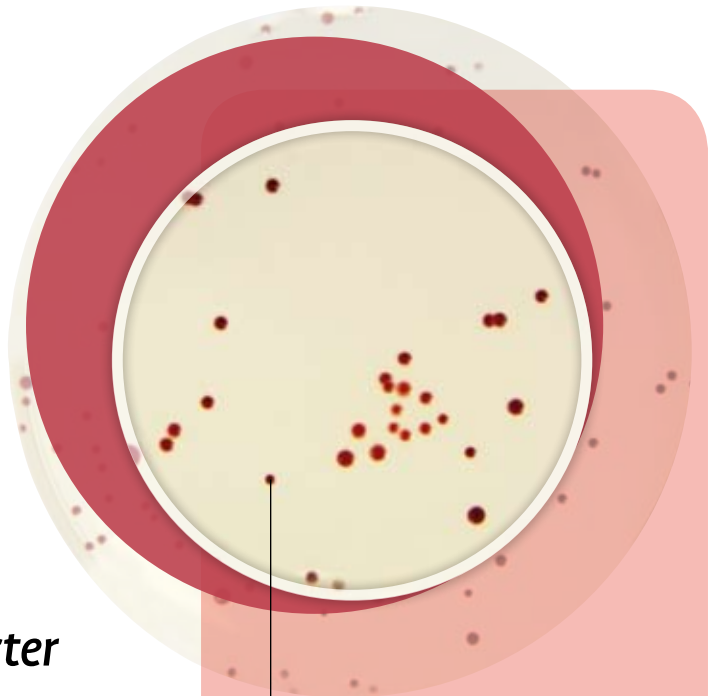

*Brick red colonie*



# CASA®

## Enumeration protocol

- Enrichment preparation (buffered peptone water or any other adapted broth)
- Spread on CASA® medium
- Incubation 44 ± 4h at 41,5°C
- Typical colonies on CASA® are red brick colonies
- Carry out confirmation test with latex agglutination test



## References

AEB140062: Bolton broth, 500g  
AEB184003/10: Bolton supplement, 10 x qty sufficient for 500 ml

AEB520420: mCCDA medium, 20 plates 90 mm  
AEB520270: CASA® medium, 20 plates 90 mm  
MGNF46: *Campylobacter* spp. latex test

### > Defrinated Horse blood

AEB300025: 25 ml  
AEB300050: 50 ml  
AEB300100: 100 ml

### > Gaspack EZ Campy. for jars

BBL60680: 20 units

### > Jars

BBL60671: Gaspack jar 18 plates - 4 liters  
BBL60672: Gaspack jar 33 plates - 6,5 liters  
BBL60673: rack jot 18 plates jars  
BBL60674: rack for 33 plates jar

### > Gaspack EZ Campy pouch system

BBL60685: 20 units



[contact@aeschemunex.com](mailto:contact@aeschemunex.com) • [www.aeschemunex.com](http://www.aeschemunex.com)

AES CHEMUNEX France  
(Headquarters)  
Rue Maryse Bastié  
Ker Lann CS17219  
F-35172 Bruz cedex  
Tel.: +33 (0) 2 2350 1212  
Fax: +33 (0) 2 2350 1200

AES CHEMUNEX, INC USA  
1 Corporate Drive  
Cranbury NJ 08512  
Tel.: +1 609 235 9272  
Fax: +1 609 235 9298

AES CHEMUNEX UK  
Endeavour House  
3<sup>rd</sup> floor • London  
Stansted Airport  
Essex • CM24 1SJ  
Tel.: +44 (0)1279 666 356  
Fax: +44 (0)1279 666 357

AES CHEMUNEX Germany  
Wilhelmstraße 49  
D-74918 Angelbachtal  
Tel.: +49(0)7265 9179 10  
Fax: +49(0)7265 9179 15

AES CHEMUNEX Spain  
Pol. Ind. Santa Margarida II  
C/ Albert Einstein, 44  
08223 TERRASSA  
(Barcelona)  
Tel.: +34 93 785 36 27  
Fax: +34 93 784 72 63

AES CHEMUNEX Italy  
Via Panà, 56/b  
35027 NOVENTA  
PADOVANA (PD)  
Tel.: +39 049 87 05 804  
Fax: +39 049 87 09 691