Twin ***sensor*** **BT**

A Rapid Test for the detection of β -Lactams & Tetracyclines in Milk.
For product code : KIT020 / KIT034 – INCUBATION TEMPERATURE : 40°C

(GB)	Instructions for use	5
(FR)	Notice d'utilisation	7
(NL)	Instructies voor gebruik	9
(D)	Gebrauchsanweisung	12
(S)	Instrucciones de uso	15
(I)	Istruzioni per l'uso	18
(TR)	Kullanma Talimatı	20
(CN)	使用说明	23
(RUS)	Инструкция по применению	24
(GR)	εγχειρίδιο χρήστη	27


Twinsensor_KIT020 / KIT034_Kit Insert_V1.0

unisensor 

Fig. A : Limits of detection

	Compounds (MRL-ppb)	Limits of detection ppb ($\mu\text{g/l}$)	Limits of detection ppb ($\mu\text{g/l}$)
		KIT020	KIT034
Penicillins	Penicillin-G (4)	2 – 3	2 – 3
	Ampicillin (4)	3 – 5	3 – 5
	Amoxicillin (4)	3 – 5	3 – 5
	Oxacillin (30)	12 – 18	12 – 18
	Cloxacillin (30)	6 – 8	6 – 8
	Dicloxacillin (30)	6 – 8	6 – 8
	Nafcillin (30)	30 – 50	30 – 50
Cefalosporins	Cefacetriple (125)	30 – 40	30 – 40
	Cefalexin (100)	> 750	> 750
	Cefalonium (20)	3 – 5	3 – 5
	Cefazolin (50)	18 – 22	18 – 22
	Cefoperazone (50)	3 – 4	3 – 4
	Cefquinome (20)	20 – 30	20 – 30
	Ceftiofur (100)	10 – 15	10 – 15
	Cephapirin (60)	6 – 8	6 – 8
Tetracyclines	Chlortetracycline (100)	30 – 40	5 – 7
	Doxycycline (100)	10 – 15	2 – 3
	Oxytetracycline (100)	50 – 60	7 – 9
	Tetracycline (100)	80 – 100	12 – 15

Fig. B : Visual Interpretation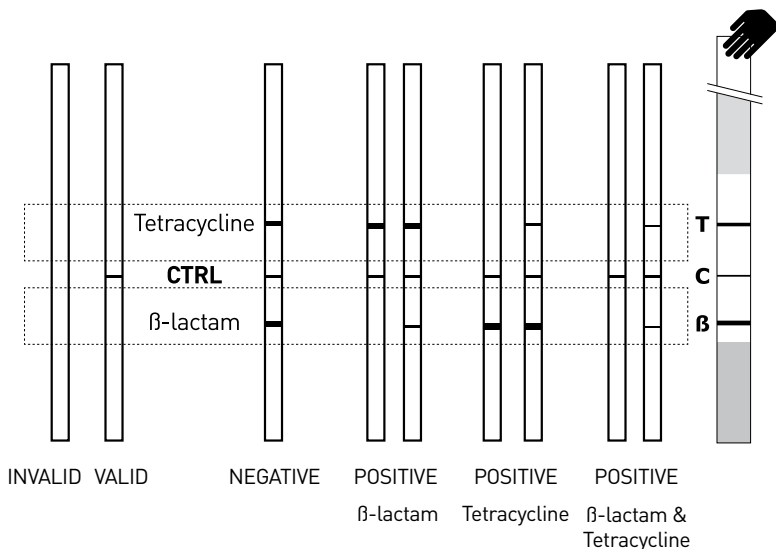


Remember this simple key :

- **T**op line : Tetracyclines
- **C**entre line : Control
- **B**ottom line : Beta Lactams

Fig. C : Positive Standard

	KIT020	KIT034
Penicillin-G	4 ppb	4 ppb
Oxytetracycline	100 ppb	20 ppb

Fig. D : How to correctly use the micropipette

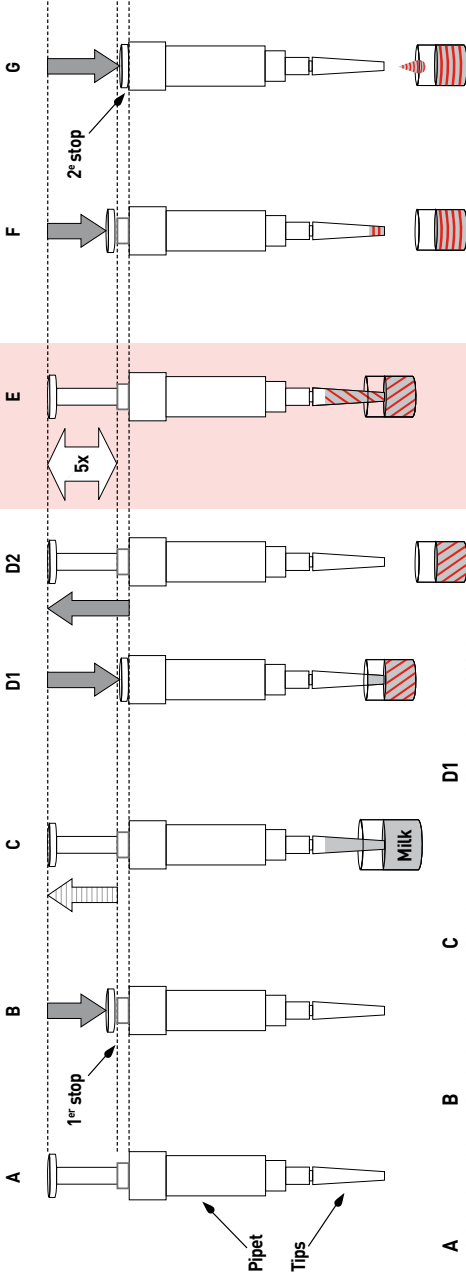- A** Original position of the pipet.
- B** Push the plunger to the 1st stop.
- C** Dip the end of the tip into the milk and slowly release allowing the plunger to return to its original position.
- D1** Transfer the 200 μ l of milk into the nominated reagent microwell and while pushing slowly the pipet plunger down to the 2nd position, pull the pipet out of the milk.
- D2** Release allowing the plunger to return to its original position.
- E** With the same tip on the pipet, mix rapidly the reagents and the milk by sucking up and rejecting the liquid 5 times consecutively so as to get a homogeneous solution.
- F&G** Eject the 200 μ l of mix by pushing the plunger to the 2nd position.

Twinsensor^{BT} code KIT020 / KIT034 are rapid tests that allow you to simultaneously detect the presence of both β -lactam and tetracycline molecules in a milk sample. **(fig. A)**

Summary of the protocol

- Add 200 μ l of milk into one reagent microwell and mix to homogeneity ;
- Incubate 3 min at 40°C ;
- Dip one Dipstick into each microwell ;
- Continue incubating for 3 min at 40°C ;
- Read the colour intensities.

Reaction mechanism

Twinsensor^{BT} is a competitive test involving two receptors in one single operation. The test requires the use of two components. The first component is a microwell containing predetermined amounts of both receptors and antibodies linked to gold particles. The second is a dipstick made up of a set of membranes with specific capture lines **(fig. B)**. For a valid test, the red control line has to be visible after the second incubation. The other two are the specific "test" lines placed on both sides of the control line. The line for β -lactam antibiotics (penicillins and cephalosporins) is located below the "control" while the line relating to tetracyclines is located above it. When the reagent from the microwell is re-suspended with a milk sample, both receptors will bind the corresponding analytes if present during the first 3-minute incubation at 40°C. Afterwards, when the dipstick is dipped into the milk, the liquid starts running vertically on the dipstick and passes through capture zones. When the sample is free of antibiotics, a colour development occurs at the specific capture lines, indicating the absence of the targeted analytes in the milk sample. On the contrary, the presence of antibiotics in the sample will not cause the coloured signal to appear at the specific capture lines.

Composition of kits

Kits coded « KIT020 / KIT034 » contain everything needed to perform 96 measurements.

- 12 pots each with 1 strip of 8 reagent microwells and 8 dipsticks ;
- 1 micropipette of 200 μ l and disposable tips ;
- « Positive Standard » containing powder to reconstitute "positive standard" raw milk of PenG and Oxytetracycline **(fig. C)**. For ease of recognition, the Positive Std is stained by a dye to give it a very pale red colour ;
- « Negative Standard » containing powder to reconstitute "negative standard" raw milk. The Negative Std is stained by a dye to give it a very pale green colour ;
- One instruction sheet.

General remarks

1. At reception, store the kit in a dry place and at a low temperature between 2°C and 8°C. Before opening, let the plastic pots reach room temperature and avoid exposure of the product to moisture and light ;
2. The standards are presented in two formats either Glass Vials or Individual Plastic Tubes. The Standard Glass vials have to be re-hydrated with 1 ml (5 X 200 μ l) of Pure Water and the Standard Individual Plastic tubes have to be re-hydrated each with 400 μ l (2 X 200 μ l) of pure Water. Mix vigorously (vortex) to avoid clots. The reconstituted glass vial can be stored in the freezer at -20°C. Do not freeze/thaw more than once ;
3. Avoid using clotted milk with Twinsensor^{BT} ;
4. The best temperature to perform the test is 40°C \pm 3°C. Use the "Heatsensor" (or alternatively a water bath). Any other type of incubator is not appropriate to perform the Twinsensor assay. (Refer to the used Heatsensor manual for setting at right temperature and timing) ;
5. After the second incubation, read the result within a 15-minute time-frame. Do not attempt to interpret the result after 15 minutes ;
6. Instrumental readings must be done with Readsensar (see user manual) ;
7. When drying, the colour intensities of the lines will become sharper ;
8. When a positive result is recorded, the test result should be confirmed ;
9. Empty one pot before opening another pot.

Milk powder dilution

In an appropriate bottle, mix 10 g of milk powder with 90 ml of warm (40°C) and distilled water. For an optimal dilution, shake vigorously.

Directions for use

This procedure is described to easily run one single sample or a set of many samples. In that case, try to perform the test in cascade and avoid any delays when mixing reagent and milk but also when adding and removing dipsticks. Make sure you have the same incubation time and temperature for each sample. You shouldn't test more than 8 samples at a time and, if there are more than 3 samples, you should use a multipipette. With more than 8 samples we recommend to share series of maximum 8 samples.

1. Choose a clean and dry place to perform the test and wash and dry your hands before starting ;
2. Connect the Heatsensor (see user manual) and wait until the temperature has stabilised at 40°C. *It takes about 10 minutes for the temperature to stabilize at 40°C ;*
3. Before opening the reagents, take the kit out of the fridge and wait until the temperature of the reagents reaches the ambient temperature. Meanwhile, read the directions for use attentively ;
 - *There are two main components of the test which are : dipsticks and the freeze-dried reagents in microwells. Both are stored in the white plastic pots.*
4. Determine how many milk samples are to be tested and write on each tube an identification number ;
 - *The milk sample must be liquid and homogeneous. There can be neither clots nor sedimentation phases. The ideal temperature of the milk sample is between 4 and 20°C.*
5. Open one plastic pot and take out as many microwells as there are milk samples to be tested (the standard positive and negative milk included if necessary) ;
 - *To open a pot of dipsticks, take off the safety ring by pressing it down the pot, take off the ring and get the stopper off the pot with your thumb ;*
 - *The pot with dipsticks should always be well closed after reagents have been taken out ;*
 - *A pot with dipsticks should be emptied before another is opened ;*
 - *Be careful, if you do not intend to use all the 8 microwells, leave the set of 8 caps on the unused ones and do not tear off the strip of the eight caps but leave it on the microwells that will not be used. Do not try to separate individual caps and put them immediately back into the white pot without damaging the dipsticks, close and make sure it is tightly sealed.*
6. Place the microwell(s) in the heating block which shows 40°C ;
7. Place a new pipette tip on the micropipette and transfer 200 µl of milk into each of the microwells. **(fig. D)**
8. **Warning : when reagents and milk are in contact, the reaction begins. Mix quickly AND IMMEDIATELY start the timer by pressing the START (RUN)* button.** The 3-minute countdown starts ;
 - *During that first incubation, each of the receptors detects whether or not there are antibiotics in the milk sample. It takes 3 minutes for the reaction to be completed.*
9. During that time, open the same pot as before, take out as many dipsticks as there are analyses in progress and close the pot. Lay the dipsticks on a clean sheet and write down the number that matches the one of the milk sample ;
10. When the 3 minutes are over, i.e. after the sound-signal, press START (STOP)* again to stop the ringing tone and dip the corresponding dipstick into each of the microwells laid in the incubator ;
 - *Make sure you put into the corresponding microwell the dipstick with the same identification number as the one of the concerned milk sample.*
11. Start the timer for the second incubation by pressing the START (RUN)* button. The 3-minute countdown starts ;
12. When the 3 minutes are over, i.e. after the sound-signal, press START (STOP)* again to stop the ringing tone and take the dipsticks out of the microwells to lay them down on a sheet of paper ;
13. Put everything back into the box and close it hermetically. If you are not planning to perform any other test with Twinsensor within the same day, store it in a fridge at a temperature ranging from 2 to 8 °C.

**(Refer to the used Heatsensor manual – For DUO Heatsensor users, steps 9 to 12 are simplified)*

Visual interpretation of results : (fig. B)

Make visual readings as follows :

14. First check whether the central control line is present. If it is not, regard the analysis as invalid and do not start (or continue) any interpretation.
15. When the central control line can be seen, interpret the two test lines as follows :
16. Examine one test line at a time and compare the intensity of the line colour of the test line with the intensity of the line colour of the control line. (Start with the bottom line of β -lactams for example) ;
 - a. If the test line is **darker** in colour than the control line, the result is **NEGATIVE**, which means that, given the sensitivity of our test, the milk sample contains no antibiotics or antibiotics at a lower level than the value stated in the enclosed table ;
 - b. If the test line is as distinct as or lighter in colour than the control line, the result is **POSITIVE**, which means that, given the sensitivity of our test, the milk sample contains antibiotics at or above the detection value stated in the enclosed table.
17. When you have interpreted one test line, do the same for the other line (continue with the upper line of Tetracycline for example) ;
18. If you hesitate, regard the sample as **POSITIVE** and confirm your interpretation by performing a second visual reading within the next 15 minutes ;
19. Write down your assessment on each of the dipsticks ;
20. Dipsticks can be archived as a permanent record if required by removing the sample pad and allowing to dry before storage. N.B. Line colour intensity will darken on drying.

If you have a Readsensor, you should read the dipstick within 15 minutes of performing the test. See the user manual of Readsensor.

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